DOS K 85

en	Fitting instructions DOS module
lt	Montavimo instrukcija DOS modulio
lv	Montāžas instrukcija Dozēšanas modulis
nl	Montagehandleiding DOS-module
no	Monteringsveiledning DOS-modul
pl	Instrukcja montażu Moduł DOS
pt	Instruções de montagem Módulo DOS
ru	Инструкция по монтажу DOS-модуль
sk	Návod na montáž DOS modul
sl	Navodila za montažo Dozirni modul
sv	Monteringsanvisning DOS-modul
tr	Montaj Talimatı DOS-Modülü
zh	安装指南 DOS 模块

en 4

lt 14

lv 24

nl 34

no 44

pl 54

pt 64

ru 74

sk 87

sl 97

sv107

tr117

zh127

Guide to the appliance	4
Caring for the environment	6
Disposal of the packing material	6
Disposing of your old machine.....	6
Installation	7
Schematic diagram showing the DOS module	7
Installing the DOS module.....	7
Suspended installation:	7
Flat installation:	7
Installation in a neighbouring unit:	8
Connection	9
Connecting the DOS module to the washer-disinfector	9
Flowmeter.....	9
Strain relief for dispensing hoses	9
Connecting the dispensing hoses.....	10
Colour coding on the dispensing hoses.....	11
Maintenance.....	11
Technical data	12

en - Guide to the appliance

The DOS module is an external dispensing system for connection to a Miele washer-disinfector only. It automatically dispenses liquid agent such as cleaning and disinfecting agent, neutralising agent and rinsing agent at the appropriate stage of the programme. This eliminates, for instance, the need to refill the dispenser in the door of your washer-disinfector. Depending on machine, up to two external dispensing systems can be connected to your washer-disinfector.

DOS modules can be retrofitted and connected if required.

 The DOS module K 85 must only be used for applications described in these instructions. Any other use, conversion or alteration is not permitted and potentially dangerous.

Miele cannot be held liable for damage resulting from incorrect or improper use or operation.

The DOS module has to be commissioned before first use and users trained in its use by a Miele technician or authorised person only.

Please observe the operating instructions for your washer-disinfector and pay particular attention to the Warning and Safety instructions regarding the use and handling of chemical agents.

Please read the installation instructions carefully before installing the DOS module. This is to ensure your safety and to prevent the risk of damage to the DOS module.


Keep these instructions in a safe place for reference.

Areas of application

DOS module K 85 is suitable for use with the following washer-disinfectors:

- PG 8562
- PG 8581
- PG 8582
- PG 8583
- PG 8591
- PG 8592
- PG 8593

Symbols on the DOS module

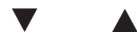Liquid agent flow direction

Warning:
Danger of electric shock


Disposal of the packing material

The packaging is designed to protect against transportation damage. The packaging materials used are selected from materials which are environmentally friendly for disposal and should be recycled.

Recycling the packaging reduces the use of raw materials in the manufacturing process and also reduces the amount of waste in landfill sites.

Disposing of your old machine


Electrical and electronic machines often contain valuable materials. They also contain materials which, if handled or disposed of incorrectly, could be potentially hazardous to human health and to the environment. They are, however, essential for the correct functioning of your machine. Please do not therefore dispose of it with your household waste.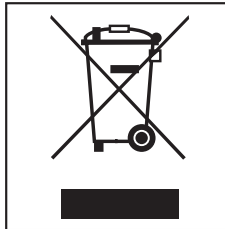

Please dispose of it at your local community waste collection / recycling centre or contact your dealer for advice.

Ensure that it presents no danger to children while being stored for disposal.


Schematic diagram showing the DOS module

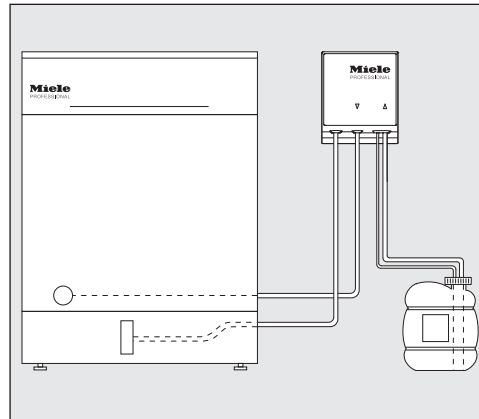Installing the DOS module

The DOS module is supplied with two screws (4.2 x 25 mm) and two plugs (6 mm) for installation.

- Select a suitably accessible position, e.g. on an adjacent wall or in a unit for installing the DOS module.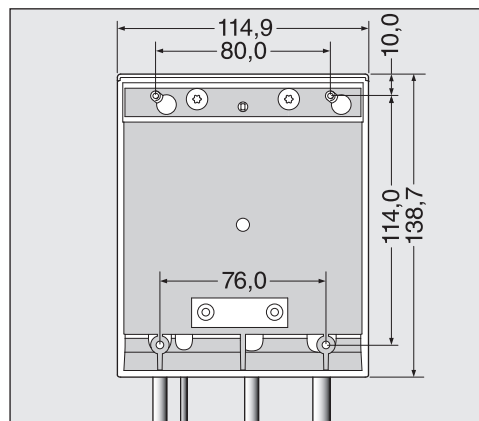

Suspended installation:

- Using the drawing as a guide fit the screws the correct distance apart (80.0 mm) and suspend the DOS module from these.

Flat installation:

Two slots, 76.0 mm apart are provided at the bottom of the DOS module for fixing the DOS module to a thin carrier or plate. 3.9 mm diameter screws (not supplied) are required for securing the DOS module.

- Drill two holes through the plate and screw the DOS module firmly in place from underneath.


en - Installation

Installation in a neighbouring unit: If more than one DOS module is being used, or if the liquid agent containers are to be stored with the DOS module, a suitable unit is available from Miele.

Make sure hoses are laid without kinks.

Connecting the DOS module to the washer-disinfector


Flowmeter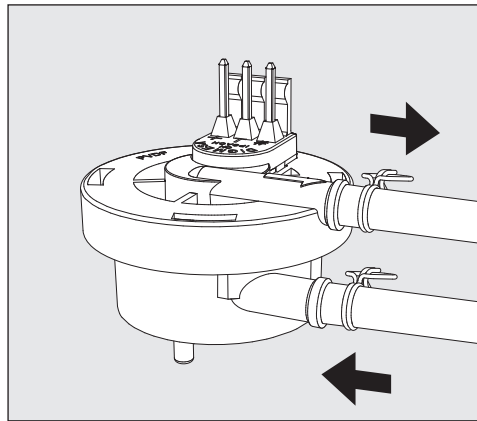

DOS module K 85 is supplied with a flowmeter which monitors the dispensing of liquid media.

Highly viscous chemical agents can cause problems with dispenser monitoring. In this case the dispenser monitoring system for laboratory applications should be switched off by Miele Service.

Strain relief for dispensing hoses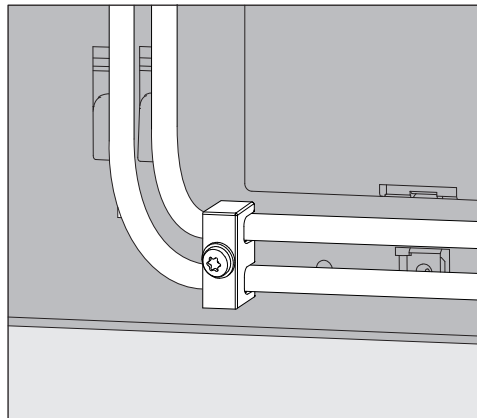

The strain relief provided with the DOS module for the dispensing hoses is fitted to the bottom right hand side of the back of the washer-disinfector to prevent the hoses from dislodging themselves from the dispenser connections during use.

The flowmeter and strain relief must only be fitted and commissioned by a Miele technician or approved person.

en - Connection


Connecting the dispensing hoses

The DOS module is tested at the factory for functionality. Any residues of the harmless test medium are a result of this testing and not evidence of previous use.

- Pull the safety cap off the end of the hose and remove the protective foil from the siphon.
- Connect the hose to the dispenser connection on the washer-disinfector. The flowmeter has to be fitted between the DOS module and the dispenser connection.
- Connect the module to the electricity supply.

Connections for the electricity supply and the dispensing hose are located at the back of the washer-disinfector. (These are described in detail in the Operating instructions for your washer-disinfector).

- Fill or replace the container, fit the siphon and secure the lid.

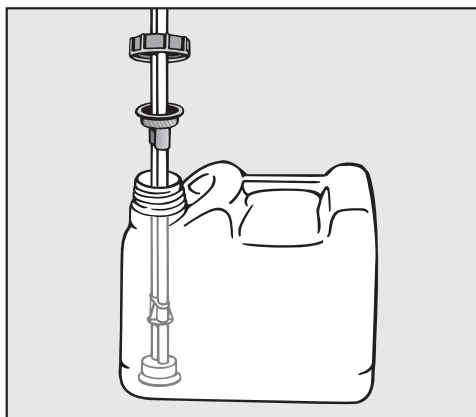- Feed the siphon into the container until it reaches the bottom.

Instructions for venting the dispensing system and setting the dosage are given in the Operating instructions for your washer-disinfector.

Colour coding on the dispensing hoses

To differentiate between the containers when using more than one DOS module, the hoses must be colour coded using the coloured labels supplied.

- Stick the label near the siphon and one of the same colour close to the free end of the hose to enable clear identification.

Miele use and recommend the following for DOS K 85:

- blue: for cleaning agent
- red: for neutralising agent
- green: for disinfecting agent / additional second cleaning agent
- white: for acidic chemical agent
- yellow: for special agents in laboratory applications

Maintenance

The dispensing system should be serviced at the same time as the washer-disinfector by Miele Service.

en - Technical data

Height	138.7 mm	
Width	114.9 mm	
Depth	114.2 mm	
Weight (net)	1.5 kg (incl. siphon, hose and cable)	
Connection	Connection to washer-disinfector	
Voltage, Connected load, Fuse rating	200 - 230 V 50 Hz 12 W 16 A	200 - 240 V 60 Hz 10 W 16 A
Throughput (nominal)	55 ml/min	
Ambient temperature	5 °C to 40 °C	
Relative humidity maximum linear decreasing to	80 % for temperatures up to 31 °C 50 % for temperatures up to 40 °C	
Installation height	as per washer-disinfector	
Pollution degree	2	
Safety class (according to DIN EN 60529)	IP21	
Test certificates awarded	EMC, VDE, radio and television suppressed	
CE-mark	2014/35/EC Low voltage directive	
Manufacturer's address	Miele & Cie. KG, Carl-Miele-Straße 29, 33332 Gütersloh, Germany	

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