---

## DOS K 85 flex

<b>en</b>	Fitting instructions Dispensing module
<b>it</b>	Montavimo instrukcija DOS modulis
<b>lv</b>	Montāžas instrukcija DOS modulis
<b>nl</b>	Montagehandleiding DOS-module
<b>no</b>	Monteringsveiledning DOS-modul
<b>pl</b>	Instrukcja montażu Moduł DOS
<b>pt</b>	Instruções de montagem Módulo DOS
<b>ru</b>	Инструкция по монтажу DOS-модуль
<b>sk</b>	Návod na montáž DOS modul
<b>sl</b>	Navodila za montažo Dozirni modul
<b>sv</b>	Monteringsanvisning DOS-modul
<b>tr</b>	Montaj Talimatı DOS Modülü
<b>zh</b>	安装说明 配给模块

<b>en</b> .....	4
<b>lt</b> .....	13
<b>lv</b> .....	22
<b>nl</b> .....	31
<b>no</b> .....	40
<b>pl</b> .....	49
<b>pt</b> .....	58
<b>ru</b> .....	67
<b>sk</b> .....	80
<b>sl</b> .....	89
<b>sv</b> .....	98
<b>tr</b> .....	107
<b>zh</b> .....	116

---

<b>Product description</b> .....	4
<b>Caring for the environment</b> .....	6
Disposal of the packing material .....	6
Disposing of your old machine.....	6
<b>Installation</b> .....	7
Schematic diagram showing the dispensing module .....	7
Fitting the dispensing module .....	7
Wall mounted: .....	7
Surface mounted:.....	7
Installation inside an adjacent unit: .....	8
<b>Connection</b> .....	9
Connecting the dispensing module to the cleaning machine .....	9
Flowmeter.....	9
Strain relief for dispensing hoses .....	9
Connecting the dispensing hoses.....	9
Colour coding on the dispensing hoses.....	10
Maintenance.....	10
<b>Technical data</b> .....	11

## en - Product description

---

The dispensing module is an external dispensing system that is intended solely for connection to Miele cleaning machines. It dispenses liquid process chemicals (cleaning agent, disinfection solution, neutralising agent or rinsing agent, for example) automatically at the appropriate stage of a programme. The dispensing module does away with the need to fill the dispenser containers in the door of the machine (if there are any). Up to two external dispensing systems can be connected depending on the type of machine.

Suitable process chemicals are available from Miele Professional.

Dispensing modules can be retrofitted and connected if required.

 Risk posed by conversions or modifications

Please note that the K 85 dispensing module must only be used for applications described in these instructions.

Any other use, conversion or modification is not permitted and potentially dangerous.

Miele cannot be held liable for damage resulting from incorrect or improper use or operation.

Initial commissioning will be undertaken by the Miele Customer Service Department or authorised service technicians, who will provide you with operating instructions.

Please follow the operating instructions for your cleaning machine and pay particular attention to warning and safety instructions when using and handling chemical agents.

 Damage caused by noncompliance with the fitting instructions.

Please read these fitting instructions carefully before fitting the dispensing module.

This prevents both personal injury and damage to the dispensing module.


Keep these fitting instructions in a safe place.

### Areas of application

DOS module K 85 is suitable for use with the following cleaning machines:

- PG 8562
- PG 8581
- PG 8582
- PG 8583
- PG 8591
- PG 8592
- PG 8593

### Symbols on the DOS module

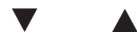Liquid agent flow direction

Warning:  
Danger of electric shock


### **Disposal of the packing material**

The packaging is designed to protect against transportation damage. The packaging materials used are selected from materials which are environmentally friendly for disposal and should be recycled.

Recycling the packaging reduces the use of raw materials in the manufacturing process and also reduces the amount of waste in landfill sites.

### **Disposing of your old machine**


Electrical and electronic machines often contain valuable materials. They also contain materials which, if handled or disposed of incorrectly, could be potentially hazardous to human health and to the environment. They are, however, essential for the correct functioning of your machine. Please do not therefore dispose of it with your household waste.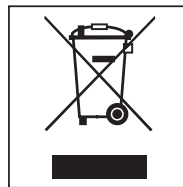

Please dispose of it at your local community waste collection / recycling centre or contact your dealer for advice.

Ensure that it presents no danger to children while being stored for disposal.

## Schematic diagram showing the dispensing module

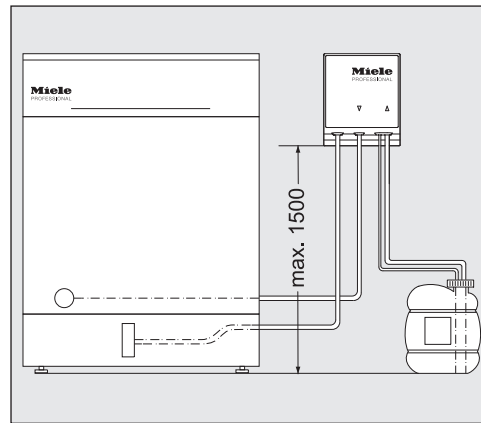
⚠ Dispensing problems caused by fitting the canister incorrectly. The canister must not be placed on top of or above the machine. Make sure that the dispensing hose is not kinked or trapped. Place the canister on the floor next to the machine or inside an adjacent unit.

## Fitting the dispensing module

The dispensing module is supplied with two screws (4.2 x 25 mm) and two plugs (6 mm) for fitting.

- Select a suitably accessible position, e.g. on an adjacent wall or inside a unit, for fitting the dispensing module.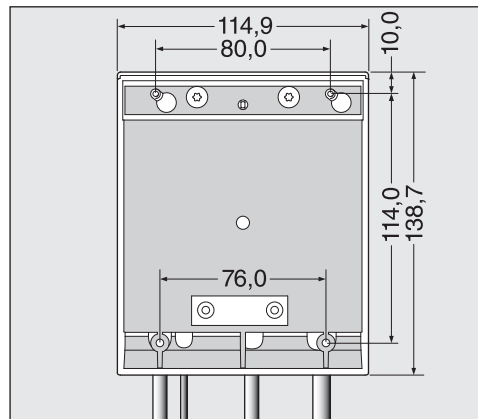

**Wall mounted:**

- Using the drawing as a guide, fit the screws the correct distance apart (80.0 mm) and suspend the dispensing module from these.

**Surface mounted:**

Two slots 76.0 mm apart are provided at the bottom of the dispensing module for fixing the module to a thin carrier or plate. Screws 3.9 mm in diameter (not supplied) are required to secure the module.

- Drill two holes through the plate and screw the module firmly in place from underneath.

## en - Installation

---


**Installation inside an adjacent unit:** If more than one dispensing module is being used, or if the liquid agent canisters are to be stored with the module, a suitable unit is available from Miele.

Make sure hoses are laid without kinks.

## Connecting the dispensing module to the cleaning machine


### Flowmeter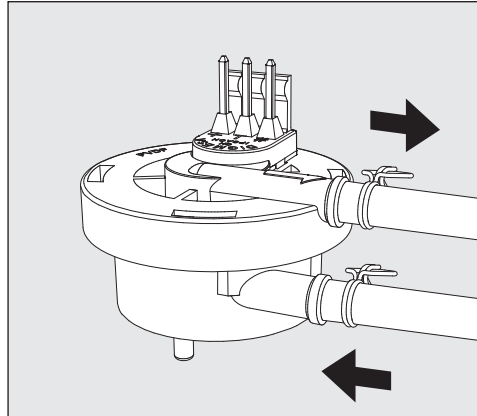

The K 85 dispensing module is supplied with a flowmeter which monitors the dispensing of liquid media.

Highly viscous chemical agents can cause problems with dispenser monitoring. In this case the dispenser monitoring system for laboratory applications should be switched off by the Miele Customer Service Department.

### Strain relief for dispensing hoses

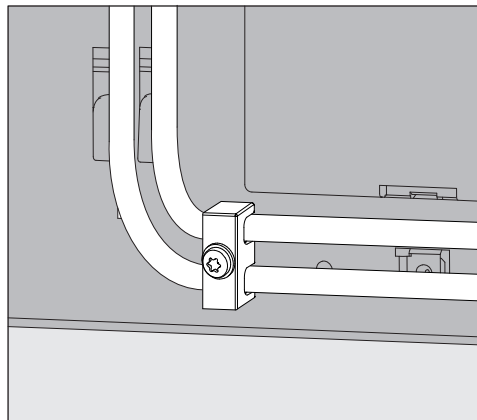The strain relief provided with the dispensing modules for the dispensing hoses is fitted to the bottom right-hand side of the back of the machine to prevent the hoses from dislodging themselves from the dispenser connections during use.

### Connecting the dispensing hoses


The dispensing module is tested at the factory for functionality. Any residues of the harmless test medium are a result of this testing and not evidence of previous use.

- Pull the safety cap off the end of the hose and remove the protective foil from the siphon.
- Connect the hose with the dispenser connection to the cleaning machine.

## en - Connection

- Connect the module to the machine's power supply.

Connections for the dispensing hoses are located on the back of the cleaning machine. These are described in detail in the operating instructions for your cleaning machine.

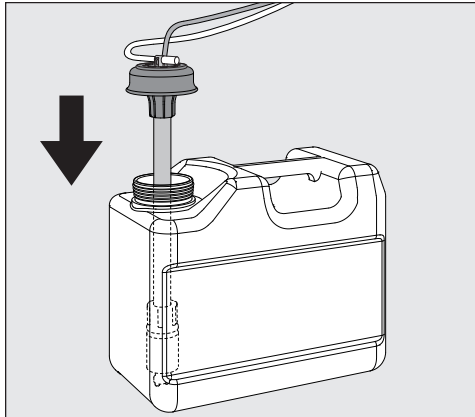- Fill or replace the container, fit the siphon and secure the lid.
- Feed the siphon into the canister until it reaches the bottom.

Follow the instructions for venting the dispensing system and setting the dosage in the operating instructions.

### Colour coding on the dispensing hoses

To differentiate between the containers when using more than one DOS module, the hoses must be colour coded using the coloured labels supplied.

- Stick the label near the siphon and one of the same colour close to the free end of the hose to enable clear identification.

Miele use and recommend the following for DOS K 85:

- blue: for cleaning agent
- red: for neutralising agent
- green: for disinfecting agent / additional second cleaning agent
- white: for acidic chemical agent
- yellow: for special agents in laboratory applications

### Maintenance

The dispensing system should be serviced at the same time as the cleaning machine by Miele Service.

## en - Technical data

Height	138.7 mm	
Width	114.9 mm	
Depth	114.2 mm	
Weight (net)	1.5 kg (incl. siphon, hose and cable)	
Connection	Connection to cleaning machine	
Voltage, Connected load, Fuse rating	200 - 230 V 50 Hz 12 W 16 A	200 - 240 V 60 Hz 10 W 16 A
Throughput (nominal)	55 ml/min	
Ambient temperature	5 °C to 40 °C	
Relative humidity maximum linear decreasing to	80 % for temperatures up to 31 °C 50 % for temperatures up to 40 °C	
Installation height	as per cleaning machine	
Pollution degree	2	
Safety class (according to DIN EN 60529)	IP21	
Test certificates awarded	EMC, VDE, radio and television suppressed	
CE-mark	2014/35/EC Low voltage directive	
Manufacturer's address	Miele & Cie. KG, Carl-Miele-Straße 29, 33332 Gütersloh, Germany	

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