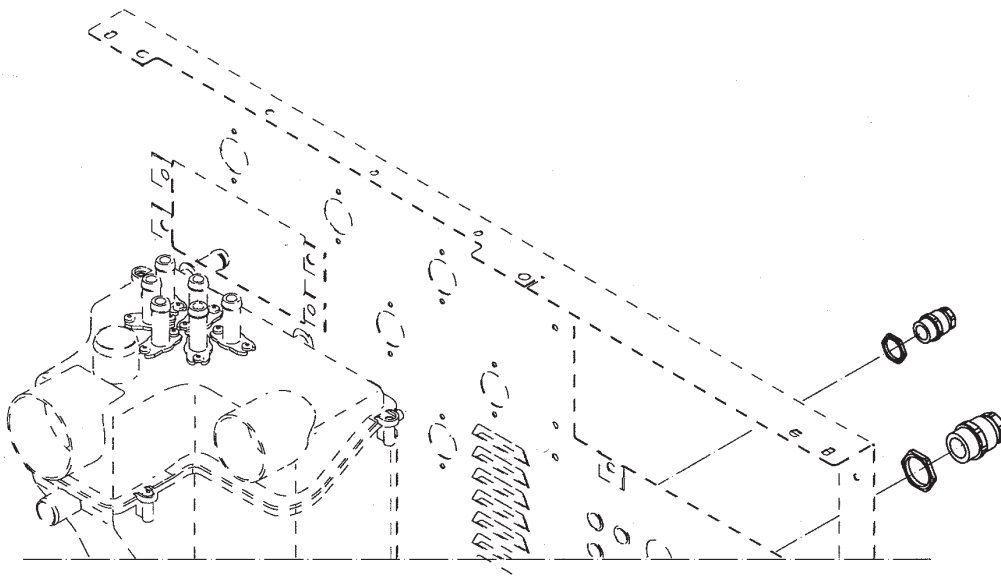
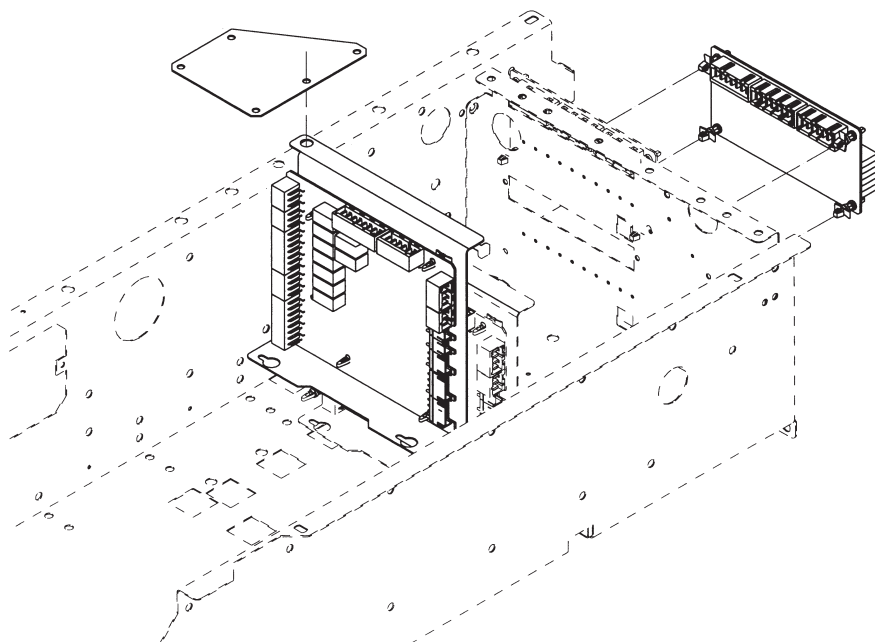
PW 6241, PW6321

<b>de</b>	Montageanweisung Modul 2 Flüssigdosierung 7-12	<b>en</b>	Fitting instructions - Module 2, liquid dispensing 7 - 12	<b>fr</b>	Notice de montage module 2 dosage liquide 7-12
<b>ni</b>	Montage-instructie moduul 2 vloeibarezeepdosering 7-12	<b>da</b>	Monteringsanvisning modul 2 dosering af flydende middel 7-12	<b>no</b>	Monteringsveiledning - Modul 2 for flytende dosering 7-12
<b>sv</b>	Monteringsanvisning: modul 2 dosering av flytande medel 7-12	<b>fi</b>	Asennusohje – Moduuli 2 nestemäisten aineiden annostus 7-12	<b>it</b>	Istruzioni di montaggio modulo 2 dosaggio liquido 7-12
<b>es</b>	Instrucciones de montaje módulo 2 Dosificación líquida 7-12	<b>am</b>	Installing Module 2 for Liquid Dispensing 7-12		

**1**

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**Note**

This **service and repair** work should only be carried out by a suitably qualified electrician (with specialist training, knowledge and experience, and recent related work experience) in accordance with all appropriate local and national safety regulations.

**Commissioning, servicing, modification, testing and maintenance** of electrical appliances should only be carried out in accordance with all appropriate legal requirements, accident prevention regulations and valid standards. All regulations of the appropriate utility supply companies and standards relating to safety (not limited to electrical safety) are to be complied with.

**Danger!**

Even with the machine switched off, mains voltage may be applied to some components.

Before any service work is commenced, the machine must be disconnected from the mains. Suitable measurements must be made to ensure that this is the case.

A general visual check and risk evaluation should always be carried out.

Incorrect conversion or service work can be dangerous (risk of fire, electric shock, etc.).

**Danger!**

Correct earthing function can be deactivated if casing parts are incorrectly assembled.

Correct earthing function must be ensured when refitting the casing.

Carry out appropriate electrical safety checks.

**Danger!**

During fitting, all accessible conductive parts that are not connected to the earth must not have any link to earth (strengthened double insulation).

To confirm this, insulation resistance and touch current measurements must be carried out.

**Note**

The kit, Mat. no. 06197490 for PW 6241, PW 6321 contains the following:

- 1 electronic module EZL 230(a), module 2 (liquid dispensing 7-12)
- 1 adapter EV 230
- 1 electronic module holder
- 8 spacers (for electronic module EZL 230(a) fastening)
- 1 coding plug MFM 2-5
- 1 wiring harness 48X-41X
- 1 wiring harness 49X-46X
- 1 wiring harness 51X-52X-101X
- 1 wiring harness 47X-100X
- 3 connection wires, green-yellow 0.75 x 240
- 1 accessory module
- 1 wiring diagram module 2, Mat. no. 09021420
- 1 wiring diagram Mat. no. 09882800
- Fitting instructions - Module 2, liquid dispensing 7 - 12, Mat. no. 06572532.

**Note**

The kit, Mat. no. 09885130 for PW 6241, PW 6321 contains the following:

- 1 electronic module EZL 230(a), module 2 (liquid dispensing 7-12)
- 1 adapter EV 230
- 1 electronic module holder
- 8 spacers (for electronic module EZL 230(a) fastening)
- 1 coding plug MFM 2-5, AWG
- 1 wiring harness 48X-41X, AWG
- 1 wiring harness 49X-46X, AWG
- 1 wiring harness 51X-52X-101X, AWG 20
- 1 wiring harness 47X-100X, AWG 20
- 1 connection wire, green-yellow 240 x 96, AWG 18
- 1 accessory module
- 1 wiring diagram module 2, Mat. no. 09021420
- 1 wiring diagram Mat. no. 09882800
- Fitting instructions - Module 2, liquid dispensing 7 - 12, Mat. no. 06572532.

List of illustrations:

- Fig. 1: Fitting cable throughfeeds on connection strip
- Fig. 2: Fitting electronic module EZL 230(a)
- Disconnect the machine from the gas (if applicable) and electric mains and ensure utilities cannot be switched on again in error.
- Remove 1 screw from each side of the lid at the top of the machine.
- Lift the front of the lid slightly, slide it to the rear and thus release the centring cones at the rear of the lid from the connection strip.
- Remove the lid upwards.
- Fit the electronic module EZL 230(a), adapter EV 230, Fig. 2, and cable throughfeeds, Fig. 1.
- Connect the electronic module EZL 230(a) and adapter EV 230, Fig. 2, in accordance with the wiring diagram.
- Modify the connection for EZL 230 46X from electronic module 2N1L to 3N1L.

**Lay and connect the wiring harnesses as follows:**

**Warning!**

The power and control wiring harnesses must be laid separately in order to avoid interference in the control circuit (electromagnetic compatibility).

Excess cable should be gathered such that current flows in opposite directions and tied in order to avoid any antenna effect.

**Connect the coding plugs as follows:**

- Connect coding plug MFM 2-5 to module 2 (liquid dispensing 7-12) **EZL 230(a)** - 42X, PIN 2-5.

**Register module 2 (liquid dispensing 7 - 12) as follows:**

- Via Service level ⇒ Main menu ⇒ Service settings ⇒ Module registration ⇒ EZL with coding 2 ⇒ Module registration? ⇒ OK button ⇒ Activate the disk symbol to save the setting.
- Switch on the dispensing display.

Via Service level ⇒ Main menu ⇒ Service settings ⇒ Basic settings ⇒ Dispensing ⇒ Dispensing display: Off ⇒ On ⇒ OK button ⇒ Back button (to go back one level) ⇒ Activate the disk symbol to save the setting.

**Note**

If the dispensing display is activated, the operator can activate and deactivate dispensing. If the dispensing display is not activated, dispensing activation and deactivation is controlled by the programme in operation only.

- To quit the service mode, switch off the machine.

**Register the dispenser pump as follows:**

- If the pump capacity (flow rate) is known, enter the figure directly in the supervisor level as follows (e.g. dispenser pump 7).  
Supervisor level ⇒ Main menu ⇒ Settings ⇒ Dispensing capacity ⇒ Dispenser pump 7: No ⇒ .....ml/min ⇒ OK button ⇒ Back button (until disk symbol is shown) ⇒ Activate the disk symbol to save the setting.
- If the pump capacity (flow rate) is unknown, calibrate the dispenser pump in the supervisor level as follows (e.g. dispenser pump 7).  
Supervisor level ⇒ Main menu ⇒ Settings ⇒ Dispenser calibration ⇒ Dispenser pump 7:  
Place the dispenser pump suction tube in a container with a level mark.  
Fill the container to the level mark.  
Press button 0/1 (dispenser pump operates for 60 s).  
Use a measuring vessel to refill the container to the level mark.  
Note the dispensed quantity via the measuring vessel and enter this figure in the display (V:.....ml).  
Activate the disk symbol to save the setting.
- Register the empty signal as follows (e.g. dispenser pump 7).  
Supervisor level ⇒ Main menu ⇒ Settings ⇒ Dispenser empty signal ⇒ Dispenser pump 7: No ⇒ Yes ⇒ OK button ⇒ Back button (to go back one level) ⇒ Activate the disk symbol to save the setting.
- Check dispensing in the programmes and carry out adjustments if necessary.
- Carry out a functional check.
- To quit the service mode, switch off the machine.

**Carry out final work:**

- Refit the machine lid.
- Carry out appropriate safety checks.
- Reconnect the machine to the electrical supply and, if appropriate, the gas supply. Switch the utilities on.

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**Note**

This service and repair work should only be carried out by a suitably qualified electrician (with specialist training, knowledge and experience, and recent related work experience) in accordance with all appropriate local and national safety regulations.

Commissioning, servicing, modification, testing and maintenance of electrical appliances should only be carried out in accordance with all appropriate legal requirements, accident prevention regulations and valid standards. All regulations of the appropriate utility supply companies and standards relating to safety (not limited to electrical safety) are to be complied with.

**Danger!**

Even with the machine switched off, voltage may be applied to some components.

Before starting any service work, the machine must be disconnected from the power supply. Suitable measurements must be made to ensure that this is the case.

A general visual check and risk evaluation should always be carried out. Incorrect conversion or service work can be dangerous (risk of fire, electric shock, etc.).

**Danger!**

Correct grounding function can be deactivated if casing parts are incorrectly assembled.

Correct grounding function must be ensured when reassembling the casing.

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**Danger!**

During installation, all accessible conductive parts that are not connected to ground must not have any link to ground (strengthened double insulation).

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List of illustrations:

- Fig. 1, Installing cable feed-throughs on the connection strip
- Fig. 2, Installing the EZL 230(a) electronic
- Disconnect the machine from the electrical supply and gas supply (if applicable) and ensure that power cannot be switched on again in error.
- Remove the left and right screws from the machine lid.
- Pull the lid slightly forward, push it back and unhook it from the connection strip.
- Remove the lid.
- Install the EZL 230(a) electronic, EV 230 adapter (Fig. 2) and cable feed-throughs (Fig. 1).
- Wire the EZL 230(a) electronic and EV 230 adapter (Fig. 2) in accordance with the wiring diagram.
- Connect electronic EZL 230 46X from electronic 2N1L to 3N1L.

**Arrange and connect the wiring harnesses:****Warning!**

The load and control current wiring harnesses should be separated from one another to prevent noise in the control circuit (electromagnetic tolerance).

Extra wiring should be laid in the opposite direction to avoid antenna-like behavior.

**Install the coded plug:**

- Install coded plug MFM 2-5 on module 2 (liquid dispensing 7-12) **EZL 230(a)** – 42X, pins 2-5.

**Register module 2 liquid dispensing 7-12:**

- **Service** → **Settings** → **Register module** → **EZL with coding 2** → **Register module?** → **OK** → **Diskette symbol** to save the registration.
- Switch on dispensing display.  
**Service** → **Settings** → **Basic settings** → **Dispensing** → **Dispensing display: Off** → **On** → **OK** → **Back** (to go back one level) → **Diskette symbol** to save display activation.

**Note**

Switching on the dispensing display enables the operator to select and deselect dispensing. Without dispensing display, only information regarding whether dispensing is active and the dispensing amount is shown.

- To quit service mode, switch the machine off.

**Connect the dispensing pump:**

- If dispensing pump capacity is known, this value can be entered directly in the operator level (e.g., dispensing pump 7).  
**Operator level →Main menu →Settings →Dispensing pump capacity →Dispensing pump 7: No →.....ml/min →OK →Back**  
(until diskette symbol appears) →**Diskette symbol** to save the capacity.
- If dispensing pump capacity is not known, the dispensing pump should be calibrated in the operator level (e.g., dispensing pump 7).  
**Operator level →Main menu →Settings →Dispensing calibration →Dispensing pump 7.**  
Place the dispensing pump's suction hose in a container with a fill line.  
Fill the container to the fill line.  
Press the **0/1** button (dispensing pump will run for 60 seconds).  
Use a beaker or measuring cup to refill the container to the fill line.  
Determine the dispensed amount in the beaker/cup and enter it in the display (V:.....ml).  
Press the diskette symbol to save the dispensed amount.
- Register the refill message (e.g., dispensing pump 7).  
**Operator level →Main menu →Settings →Dispenser refill message →Dispensing pump 7: No → Yes →OK →Back** (to go back one level) →**Diskette symbol** to save the refill message.
- Check dispensing in all programs and adjust as necessary.
- Carry out a functional check.
- To quit service mode, switch the machine off.

**Finish up:**

- Re-install the machine lid.
- Carry out a safety check.
- Reconnect the machine to the electrical and gas supplies and switch it on.

**am**